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SHORT PROCESSING TIMES

AVOID EXCESSIVE HEAT GENERATION AND MOTOR OVERLOAD

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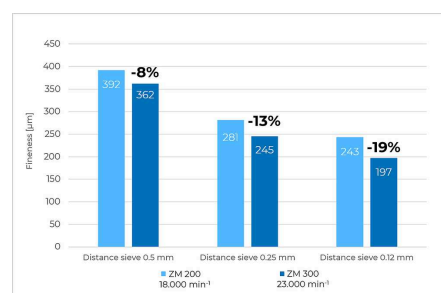
FINER END PRODUCTS AND HIGHER SAMPLE THROUGHPUT

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
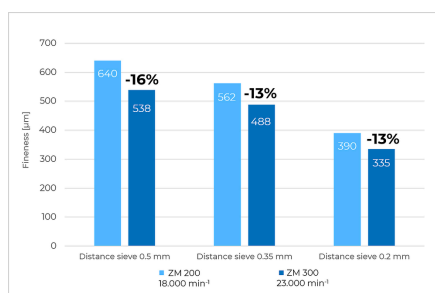
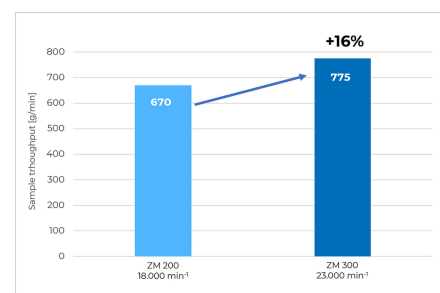
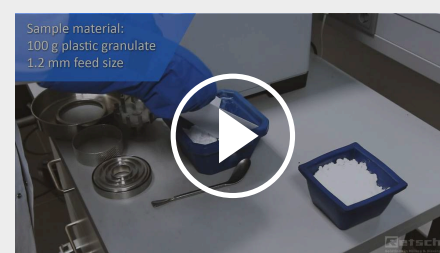
In combination with the cyclone suction unit, the Ultra Centrifugal Mill ZM 300 mill is even suitable for grinding light and heat-sensitive materials.

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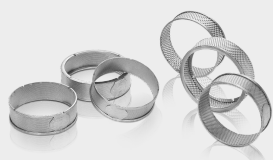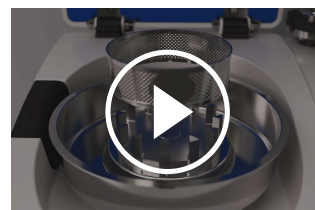
CONVENIENT OPERATION – EASY CLEANING

The large touch display with rotary knob facilitates the setting of all grinding parameters and provides current values such as cassette temperature and motor load.

Thanks to the screwless push-fit system and the patented cassette principle, grinding tools can be inserted and removed by hand. Plug-in rotors and ring sieves enable quick and easy cleaning. All components that come into contact with the product can be cleaned using running water or a dishwasher.

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The ZM 300 is suitable for grinding a wide variety of materials:

waste, graphite, collagen, drugs, fertilizers, feed pellets, grains, spices, rubber, wood, coffee beans, bones, coal, food, corn, minerals, paper, plant parts, polymers, powder coatings, rice, seeds, straw, sweets, tobacco, textiles, chemical products, electronic components, dried fruits and vegetables, dried larvae, pharmaceutical products.

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The grinding process is carried out by impact and shearing forces between the rotor and the fixed ring sieve. The sample material is fed into the rotor via a funnel with splash protection, accelerated outwards by centrifugal force, and pre-ground by the wedge-shaped rotor teeth. The subsequent fine grinding takes place in the narrow gap between the rotor and the sieve.

Thanks to this two-stage process, sample preparation is particularly gentle and fast. Since the ground material only remains in the grinding chamber for a short time, the characteristic properties of the sample are retained. The end product is collected in the collection vessel, cyclone, or paper filter bag that surrounds the grinding chamber.

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www.retsch.com/zm300

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ULTRA CENTRIFUGAL MILL ZM 300

Ultra Centrifugal Mill ZM 300 with 900 ml cassette
(please order push-fit rotor and ring sieve separately)

20.824.0001		ZM 300	230 V, 50/60 Hz
20.824.0002		ZM 300	110-120 V, 50/60 Hz

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RING SIEVES OF STAINLESS STEEL, WITH REINFORCED RIM

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ACCESSORIES ZM 300 FOR ABRASIVE PRODUCTS

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ACCESSORIES ZM 300 FOR GRINDING WITHOUT HEAVY-METAL CONTAMINATION

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22.355.0012 Cassette 900 ml with coating of titanium-niobium, complete (pan and lid with gasket)

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CYCLONES ZM 300

CYCLONE INCL. CASSETTE PAN WITH OUTLET AND SUPPORT

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22.748.0005		産業用掃除機, 230 V, 50/60 Hz (other electrical versions and spare filter bags available upon request)
-------------	---	---

ACCESSORIES FOR COLD GRINDING ZM 300

22.354.0007		Cryo kit for embrittlement of samples in liquid nitrogen or dry ice snow (incl. 2 insulated containers (1 and 4 liter), 1 spoon spatula, 1 pair of safety glasses)
-------------	--	---

ADDITIONAL ITEMS ZM 300

22.355.0018		Cassette complete, stainless steel, incl. pan 900 ml, lid with gasket
22.355.0021		Cassette complete, stainless steel, incl. pan 1200 ml, lid with gasket
22.355.0020		Cassette complete, stainless steel, incl. pan with outlet, lid with gasket
22.355.0012		Cassette 900 ml with coating of titanium-niobium, complete (pan and lid with gasket)
03.010.0058		Cassette pan 900 ml, stainless steel
03.010.0059		Cassette pan 1200 ml, stainless steel
02.010.0060		Cassette pan with outlet, stainless steel
03.010.0062		Cassette pan 900 ml, titanium-niobium coating
22.355.0015		Cassette lid, stainless steel, with gasket
22.355.0011		Cassette lid, titanium-niobium coating, with gasket
22.936.0009		パッキン カセット蓋用 5本組
02.011.0078		連続捕集用受器 5L ステンレス製
02.011.0031		連続試料受器 3L ステンレス製
02.107.0634		Lid for collecting receptacle, 5 litres
02.107.0493		Lid for collecting receptacle, 3 litres
02.186.0010		サイクロン用フィルタチューブ ナイロン製
01.706.0318		Labyrinth disc
03.114.0092		Gasket for labyrinth disc, 1 piece

03.785.0156		ホッパー
03.143.0049		Hopper insert for rebound safety
22.524.0006		ダストフィルタ枠 (フィルタ付き)
22.524.0007		Dust filter for ZM 300, 10 pcs.
99.200.0033		IQ/OQ Documentation for ZM 300

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70.938.0001		DR 100 drive unit	100V, 50 Hz
70.938.0002		DR 100 drive unit	100V, 60 Hz
72.020.0018		Feeding kit DR 100-75/40 for ZM 300, incl. push-fit feed chute 75/40 mm, length 250 mm, holder for push-fit feed chute, hopper 3.5 litres and fixture for hopper, stand with data cable	

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