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Professional Food Sample Preparation

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[Click to view video](#)

FOR LARGE FOOD SAMPLES

- | Efficient size reduction of up to 4500 ml feed quantity due to powerful 1.1 kW motor (peak power input >3 kW)
- | Pre- and fine grinding in one mill: cutting action in regular mode, grinding by impact in reverse mode, pre-grinding in interval mode
- | All parts in contact with the sample material are autoclavable
- | Perfect adaptation to application requirements by variable speed from 500 to 4000 min⁻¹ with an increment of 100 min⁻¹
- | Optional gravity lid for automatic reduction of chamber volume
- | 10 SOPs can be stored
- | Grinding jar in plastic and stainless steel available

CRYOGENIC GRINDING WITH DRY ICE

Tough or elastic materials which cannot be homogenized at room temperature, are safely and effectively homogenized in the knife mill GM 300 with the help of dry ice. For this application, a special cryo lid with a small opening to allow for possible overpressure to escape is available. Used in combination with a stainless-steel container and full metal knife, it ensures safe and efficient cold grinding.

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Operation of the laboratory knife mill GRINDOMIX GM 300 is exceptionally comfortable and safe. The grinding container and knife are easily attached without tools. When the set time has expired, the motor is automatically switched off and the container can be removed. Thus, the GRINDOMIX, unlike many commercial household mixers, allows filling and emptying the container outside the mill. The knife remains on the knife holder during the process and can be easily removed for cleaning afterwards. Thanks to this quick and easy procedure, cross contaminations caused by sample residues are successfully prevented. Another advantage of the GRINDOMIX knife mills: the grinding tools are autoclavable. The GM 300 features a convenient 1-button operation with graphic display. All parameters are digitally set and 10 Standard Operating Procedures (SOPs) can be stored.

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A range of different containers, lids and knives makes the GRINDOMIX GM 300 a truly universal device. The selection of the grinding container depends on the products to be processed. Standard plastic containers are suitable for the majority of applications. Other accessories are available for special applications.

| Gravity lid made of autoclavable plastic

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| Special lid for dry ice applications

Made of autoclaveable plastic, it allows possible overpressure to escape through a small opening.

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| Knives

4 different knives available: stainless steel, stainless steel with serrated blades, full metal blades & cylinder, with titanium-niobium coating for heavy-metal-free size reduction. All are autoclavable.

| Scraper

Facilitates the recovery of sticky samples from the container.

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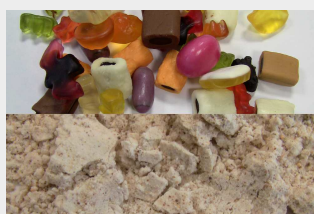
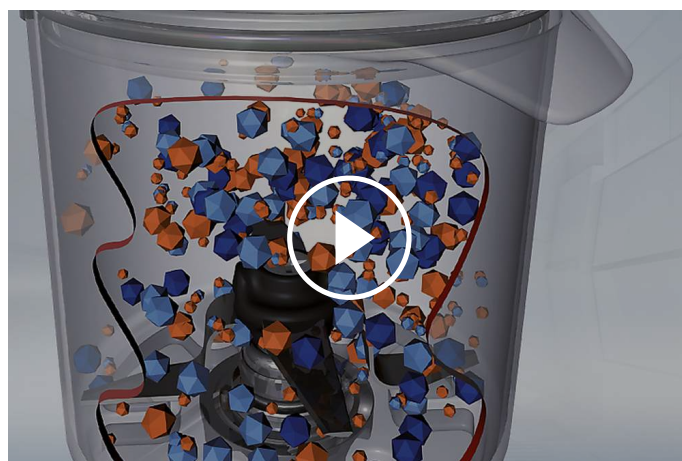
The laboratory knife mill GM 300 features four sharp, robust blades, which rotate in the center of the grinding container. Depending on the rotational direction, size reduction is effected with the blunt side (preliminary size reduction) or the sharp side (fine grinding). The knives are protected with a counterblade from damages by hard sample materials.

The knife is indirectly driven by a powerful motor of 1.1 kW which can reach temporary peaks of 3 kW. A pre-selectable speed which is maintained electronically ensures reproducible results.

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KNIFE MILL GRINDOMIX GM 300

Knife Mill GRINDOMIX GM 300 complete with grinding container 5 litre of autoclavable plastic, standard lid, knife and scraper

20.252.0001

GM 300

220-230 V, 50/60 Hz

other electrical versions available for the same price

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02.107.0522

Lid
autoclavable plastic
for dry ice applications

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02.446.0030		Knife of stainless steel, autoclavable
02.446.0054		Knife of stainless steel with serrated blades, autoclavable
02.446.0049		Knife of stainless steel (full metal blades and knife cylinder), autoclavable
02.446.0052		Knife with titanium-niobium coating, autoclavable, for grinding without heavy-metal contamination

ADDITIONAL ITEMS GM 300

05.723.0002		Съргалка
22.111.0002		Set of gaskets for lids, 10 pcs.
99.200.0011		IQ/OQ Documentation for GM 300